Inverse-square law and absorption of gamma or beta rays with the Geiger-Müller counter



Physics	Modern Physics Nuclear & particle physics		particle physics
Applied Science	Medicine	Nuclear Medicine	
Difficulty level	RR Group size	Preparation time 45+ minutes	Execution time 45+ minutes
This content can also be found online at:		×.	



http://localhost:1337/c/5fd7ea8dce3751000353f424





General information

Application

PHYWE



The understanding of the distance and absorption behaviour of radiation is very importent for the safe handeling of radioactive materials. As such this understanding is very important in nuclear power plants.



Other information (1/2)

section.

attenuation coefficient calculated.

and the counter tube.

absorbers.

PHYWE



Prior





Main

principle

The inverse square law of distance is demonstrated with the gamma radiation from a 60 Co preparation, the half-value thickness and absorption coefficient of various materials determined with the narrow beam system and the corresponding mass

The prior knowlege required for this experiment is found in the theory

Other information (2/2)

PHYWE



The goal of this experiment is to investigate the distance and absorbtion behaviour of beta and gamma radiation.

Learning

objective



Tasks

2. Determine the half-value thickness $d_{1/2}$ and the absorption coefficient μ of a number of materials by measuring the impulse counting rate as a function of the thickness of the irradiated material. Lead, iron, aluminium, concrete and Plexiglas are used as

1. Measure the impulse counting rate as a function of the distance between the source

3. Calculate the mass attenuation coefficient from the measured values.



Theory (1/5)

PHYWE

PHYWE

The cobalt isotope ${}^{60}_{27}$ Co has a half-life of 5.26 years; it undergoes beta-decay to yield the stable nickel isotope ${}^{60}_{28}$ Ni - see Fig. 1.

As with most beta emitters, disintegration leads at first to daughter nuclei in an excited state, which change to the ground state with the emission of gamma quanta. Whereas the energy levels of the beta electrons can assume any value up to the maximum because of the antineutrinos involved, the gamma quanta which participate in the same transition process have uniform energy, with the result that the gamma spectrum consists of two discrete, sharp lines (Fig. 1).



Theory (2/5)

The impulse counting rate $N^{(r)}$ per area A around a pointsource decreases in inverse proportion to the square of the distance provided the gamma quanta can spread out in straight lines and are not deflected from their track by interactions.

$$2\cdot \mathbf{r}_1 \hspace{1.5cm} A_2 = 4\cdot A_1 = \left[rac{r_2}{r_1}
ight]^2$$

The reason for this is that, as shown by Fig. 2, the area of a sphere round the source through with a beam of rays passes, increases as the square of the distance r.



Fig. 2: Law of distance relating to rays which are propagated in a straight line from a point source.

PHYWE

 $r_2 =$

www.phywe.de

PHYWE

Theory (3/5)

In vacuum (in air), therefore

$$rac{N^{(r)}}{A} = rac{N^{(0)}}{A} \cdot rac{1}{4\pi} r^{-2}$$

If we plot the counting rate $N^{(r)}$ versus the distance r on a loglog scale, we obtain a straight line of slope –2.

From the regression lines from the measured values in Fig. 3, applying the exponential expression

 $N_{(r)} = a \cdot r^b$ we obtain the value $b = -2.07 \pm \, 0.01$ for the exponent.

This thus proves the applicability of the inverse square law. The attenuation of the gamma rays when they pass through an absorber of thickness d is expressed by the exponential law.

Theory (4/5)

 $N_{(d)} = N_{(0)} \cdot e^{-\mu d}$

where N(d) is the impulse counting rate after absorption in the absorber, and N(0) is the impulse counting rate when no absorption takes place: μ is the absorption coefficient of the absorber material and depends on the energy of the gamma quantum. Fig. 3: Absorption of gamma rays by leads as a function of the energy (μ_{Co} fraction due to Compton effect, μ_{Ph} fraction due to photoelectric effect, μ_{Pa} fraction due to pair formation). The total absorption coefficient (attenuation coefficient) is $\mu = \mu_{Co} + \mu_{Ph} + \mu_{Pa}$





PHYWE

Theory (5/5)

PHYWE

The absorption of the gamma rays is brought about by three independent effects – the Compton effect, the photelectric effect and pairformation. The relative contributions of these three effects to total absorption depends primarily on the energy of the quanta and on the atomic number of the absorber (Fig. 3).

We can see from the $\frac{\mu}{E}$ curves that lead is particularly suitable as an absorber of gamma rays of low or high energy. The attenuation of gamma rays therefore takes place predominantly in the electron shell of the absorber atoms. The absorption coefficient μ should therefore be proportional to the number of electrons in the shell per unit volume, or approximately proportional to the density ρ of the material. The mass attenuation coefficient $\frac{\mu}{a}$ is therefore roughly the same for the different materials.

The half-value thickness $d_{1/2}$ of a material is defined as the thickness at which the impulse counting rate is reduced by half, and can be calculated from the absorption coefficient in accordance with

$$d_{1/2} = rac{\ln 2}{\mu}$$

PHYWE

Equipment

Position	Material	Item No.	Quantity
1	PHYWE Geiger-Müller Counter	13609-99	1
2	Geiger-Mueller counter tube, type B, BNC cable 50 cm	09005-00	1
3	Vernier calliper stainless steel 0-160 mm, 1/20	03010-00	1
4	Absorption plates for beta rays	09024-00	1
5	Base plate for radioactivity	09200-00	1
6	Counter tube holder on fixating magnet	09201-00	1
7	Source holder on fixing magnet	09202-00	1
8	Plate holder on fix. magnet	09204-00	1
9	Absorption material, lead	09029-01	1
10	Absorption material, iron	09029-02	1
11	Absorption material, aluminium	09029-03	1
12	Absorption material, concrete	09029-05	1
13	Radioactive sources, set	09047-40	1



PHYWE



Setup and Procedure

Setup

According to Fig. 4

The distance between the front edge of the source rod and the counting tube window is approximately 4 cm; consequently, the absorption plates can be easily inserted into the radiation path.





PHYWE



Evaluation

Evaluation (1/2)

PHYWE

From the regression lines from the measured values we obtain the following values for $\frac{\mu}{\rho}$ and for $d_{1/2}$ and μ/ρ , with the relevant standard errors, using the exponential expression

 $N = ae^{bd}$

Lead: $(\rho = 11.34 \ gcm^{-3})$ Aluminium: $(\rho = 2.96 \ gcm^{-3})$ $\mu = 0.62 \ cm^{-1}$ $s_{\mu} = 0.009 \ cm^{-1}$ $\mu = 0.15 \ cm^{-1}$ $s_{\mu} = 0.01 \ cm^{-1}$ $d_{1/2} = 1.12 \ cm^{-1}$ $s_{d_{1/2}} = 0.02 \ cm$ $d_{1/2} = 4.6 \ cm^{-1}$ $s_{d_{1/2}} = 0.3 \ cm$ $\frac{\mu}{\rho} = 0.055 \ cm^2 g^{-1}$ $s_{\frac{\mu}{\rho}} = 0.001 \ cm^2 g^{-1}$ $\frac{\mu}{\rho} = 0.056 \ cm^2 g^{-1}$ $s_{\frac{\mu}{\rho}} = 0.004 \ cm^2 g^{-1}$



Evaluation (2/2)

PHYWE

$$\begin{aligned} \text{Iron:} (\rho = 7.86 \ gcm^{-3}) & \text{Concrete:} (\rho = 2.35 \ gcm^{-3}) \\ \mu = 0.394 \ cm^{-1} & s_{\mu} = 0.006 \ cm^{-1} & \mu = 0.124 \ cm^{-1} & s_{\mu} = 0.009 \ cm^{-1} \\ d_{1/2} = 1.76 \ cm^{-1} & s_{d_{1/2}} = 0.03 \ cm & d_{1/2} = 5.6 \ cm^{-1} & s_{d_{1/2}} = 0.4 \ cm \\ \frac{\mu}{\rho} = 0.050 \ cm^2 g^{-1} & s_{\frac{\mu}{\rho}} = 0.001 \ cm^2 g^{-1} & \frac{\mu}{\rho} = 0.053 \ cm^2 g^{-1} & s_{\frac{\mu}{\rho}} = 0.004 \ cm^2 g^{-1} \\ \text{Plexiglas:} (\rho = 1.119 \ gcm^{-3}) \\ \mu = 0.078 \ cm^{-1} & s_{\mu} = 0.004 \ cm^{-1} \\ d_{1/2} = 8.9 \ cm^{-1} & s_{d_{1/2}} = 0.5 \ cm \\ \frac{\mu}{\rho} = 0.066 \ cm^2 g^{-1} & s_{\frac{\mu}{\rho}} = 0.003 \ cm^2 g^{-1} \end{aligned}$$

Comment

PHYWE

The procedure and evaluation are shown here in an exemplary experiment for γ -quanta; however, they can also be performed in an analogous manner for electrons. In the latter case, the Sr-90 source rod from the radioactive sources set (09047.50) and the absorption plate set for β -radiation (09024.00) must be used.

